

High Performance Cell-Free Wheat Germ Protein Expression System

INSTRUCTION MANUAL

WEPRO8240 (15N)/WEPRO8240G (15N)/WEPRO8240H (15N) and WEPRO8240 (13C/15N)/WEPRO8240G (13C/15N)/WEPRO8240H (13C/15N) Core Kit

These kits provide reagents to prepare stable-isotope-labeled proteins for protein NMR experiments:

Depending on the kit, reagents are provided for only ^{15}N labeling or ^{13}C and ^{15}N combined labeling

Product Number(s): CFS-C8N, CFS-C8GN, CFS-C8HN, CFS-C8CN, CFS-C8GCN, CFS-C8HCN

Version/date: Version 2.0_eng/March 2022

This Product has a shelf life of 1 year being properly stored at -80°C .

CFS products are for research use only.

All our manuals are available for download in PDF versions on our homepage at:
<https://www.cfsciences.com/eg/>. Work with electronic versions where possible to reduce paper waste.



Our products are produced under a strict quality management system offering high-quality reagents including wheat germ extracts from wheat obtained by natural farming in Japan.

Table of Contents

Important Information.....	3s
Shipment and Storage.....	3
Safety	3
For your convenience:	3
Introduction	4
Kit Contents.....	5
Protocol.....	7
Time Requirements.....	7
Transcription Reaction Using DNA Template.....	8
Translation Reaction Using RNA Prepared from DNA Template	9
Troubleshooting.....	11
Additional Information.....	12
Customer Information	12
Label License Policy.....	12
Trademarks	13
Others	13
Contact.....	13
Appendix	14

Important Information

Shipment and Storage

Our products are shipped on dry ice. Wheat germ extracts are temperature sensitive and must always be kept frozen. Store kit at -80°C right upon arrival and only thaw reagents when needed. Avoid repeated freeze/thawing cycles. Prepare aliquots of the wheat germ extract on first use if you want to keep making more expression experiments later; refer to the protocol below on how much extract is needed per reaction. Do not freeze/thaw the wheat germ extract more than three times.

Safety

This kit can be used in a regular molecular biology laboratory. We strongly advise to work under RNase-free conditions. Refer to a laboratory handbook for more information on how to work under RNase-free conditions.

Read the protocol carefully before starting the experiment.

Do not drink or eat in the laboratory, and always wear gloves and a lab coat while working in the lab.

Wash hands before and after doing an experiment. If you get any reagent(s) in your eyes or on your skin, wash eyes or skin immediately with water. Although this kit does not contain any hazardous reagents, do not take any risk.

Inform yourself about the necessary precautions for performing SDS-PAGE experiments using high voltage, and toxic chemicals in case you wish to prepare your own gels.

Safety Data Sheets (SDS) for our products can be downloaded from our homepage at:

<https://www.cfsciences.com/eg/>

Contact CellFree Sciences for further support and advice if you have any questions on the experiments described herein and materials provided with this product. Contact information is given at the end of this manual.

For your convenience

CellFree Sciences is providing short versions of our protocols ("Bench Notes"). Use these Bench Notes to setup your transcription and translation experiments at your workplace. They only contain the basic information needed for setting up the experiments. Use the checkmarks in the Bench Notes to assure that all pipetting steps have been completed correctly.

Introduction

The WEPRO®8240 Series NMR Core Kits provide you with the latest version of our high-performance wheat germ cell-free protein expression system for conducting protein labeling experiments. CellFree Sciences provides these labeling kits for NMR in two different versions: In the ^{15}N version all nitrogen atoms in all amino acids are ^{15}N -labeled; in the $^{13}\text{C}^{15}\text{N}$ version all carbon and nitrogen atoms in all amino acids are ^{13}C and ^{15}N -labeled. Refer to the appendix to this manual for a list on the labeled amino acids that had been used to prepare the 40xSUB-AMIX® SGC FOR NMR translation buffers.

This manual describes protocols to perform *in vitro* transcription and translation experiments at three different reaction scales using a bilayer reaction format for protein expression. CellFree Sciences provides WEPRO®8240 wheat germ extracts in three different versions that can be used under identical reaction conditions. Besides the regular WEPRO®8240 extract for universal protein expression, we have a dedicated “WEPRO®8240G Version” for the preparation of GST-tagged proteins and a dedicated “WEPRO®8240H Version” for the preparation of His-tagged proteins. Wheat germ extracts of the WEPRO®8240G and WEPRO®8240H versions have been precleared to remove wheat proteins from the extracts that would otherwise bind to glutathione or Ni resins during the purification step.

CellFree Sciences provides dedicated expression vectors optimized for use with our wheat germ cell-free protein expression system. We recommend using our expression vectors for template preparation. It is also possible to prepare expression templates by PCR methods for rapid expression testing and high-throughput studies. Note that using linear DNA templates from PCR reactions will commonly lead to lower protein yields as linear DNA is less stable than circular plasmid DNA.

This product does not provide any expression vector for the wheat germ cell-free expression system. We are offering a pEU vector set (CFS-PEU-v1.0) comprising several expression vectors optimized for use with our expression system. You can find more information on the vector set including sequence information and vector maps on our homepage at: <https://www.cfsciences.com/eg/products/vector>.

To support the formation of disulfide bonds, CFS offers the Disulfide Bond PLUS Expression Kit (CFS-EDX-DB), a version of the Premium PLUS Expression Kit, including a PDI and Ero1 α enzyme mix and altered translation reagents. The larger Disulfide Bond Enhancer Enzyme Set (CFS-DBEE) provides the PDI and Ero1 α enzyme mix and modified buffers. A version of the WEPRO8240H Core Kit (CFS-C8H/CFS-C8HN/CFS-C8HCN) is required for working with the Disulfide Bond Enhancer Enzyme Set.

Refer to our homepage or contact our support team for more information on how to use our cell-free protein expression system. The contact information is given at the end of the manual.

Kit Contents

The following table summarizes the reagents provided with this product:

Item	Quantity	Volume	Vial	Vial Color
WEPRO®8240/WEPRO®8240G/ WEPRO®8240H Extract	1	1000 µl	2.0 ml tube	Clear
5xTranscription Buffer LM	1	240 µl	1.5 ml tube	Yellow
NTP Mix (25 mM)	1	120 µl	1.5 ml tube	Blue
RNase Inhibitor (80 U/µl)	1	15 µl	1.5 ml tube	Orange
SP6 RNA Polymerase (80 U/µl)	1	15 µl	1.5 ml tube	Green
Creatine Kinase (20 mg/ml)	1	20 µl	1.5 ml tube	Red
40xSUB-AMIX® SGC FOR NMR S-1*	1	600 µl	1.5 ml tube	Purple
40xSUB-AMIX® SGC FOR NMR S-2*	1	600 µl	1.5 ml tube	White
40xSUB-AMIX® SGC FOR NMR S-3*	1	600 µl	1.5 ml tube	Brown
40xSUB-AMIX® SGC FOR NMR S-4*	1	600 µl	1.5 ml tube	Gray

Refer to the table below for more information on how to handle and store the reagents:

Item	Description	Storage
WEPRO®8240/ WEPRO®8240G/ WEPRO®8240H	WEPRO®8240/WEPRO®8240G/WEPRO®8240H wheat germ extracts are temperature sensitive! Immediately after thawing the extracts place them on ice. Upon thawing for the first time, prepare aliquots for later use and store them at -80°C. Do not freeze/thaw wheat germ extract more than three times. We recommend using liquid nitrogen to freeze the extracts.	-80°C
40xSUB-AMIX® SGC FOR NMR (*1) (S-1, S-2, S-3, S-4)	The 40xSUB-AMIX® SGC FOR NMR translation buffer is provided in four separate master mixes (S-1, S-2, S-3, S-4). It is possible to prepare onetime a large volume of the 2xSUB-AMIX® SGC FOR NMR translation buffer, and to store aliquots at -80°C for later use. Prepare 6 ml of nuclease-free water and add 550 µl of each master mix to the water, then adjust the final volume to 11 ml for making the 2x translation buffer. Do not mix the 40xSUB-AMIX® SGC FOR NMR translation buffers directly as some precipitate may occur. Avoid unnecessary freeze/thawing of the 2xSUB-AMIX® SGC FOR NMR translation buffer.	-80°C
5xTranscription Buffer LM	After thawing for the first time, divide 5xTranscription Buffer LM into appropriate aliquots convenient for later use. The buffer can be stored at -20°C.	-20°C
NTP Mix	The NTP Mix contains ATP, GTP, CTP, and UTP at a concentration of 25 mM. The NTP Mix can be stored at -20°C.	-20°C
SP6 RNA Polymerase	Provided in 50% glycerol, and can be stored at -20°C.	-20°C

INSTRUCTION MANUAL

RNase Inhibitor	Provided in 50% glycerol, and can be stored at -20°C.	-20°C
Creatine Kinase (*2)	Creatine Kinase is temperature sensitive, and we recommend to always use fresh enzyme preparations. The Creatine Kinase in the kit is provided at a concentration of 20 mg/ml. This concentration can be directly used to set up translation reactions on a MEDIUM and LARGE scale. The enzyme must be diluted with nuclease-free water to 1 mg/ml for setting up SMALL scale translation reactions.	-80°C

*1: Depending on the kit, the 40xSUB-AMIX® SGC FOR NMR buffers are provided as a (¹⁵N) or (¹³C/¹⁵N) versions. In this manual, both versions are referred to as “SUB-AMIX® SGC FOR NMR.” Refer to the tables in the Appendix to this manual for a list on the labeled amino acids used to prepare each buffer.

*2: Creatine Kinase can be purchased from Sigma-Aldrich, Catalog No. 10127566001. This will provide you with the same enzyme we had added to the kit; we cannot comment on the use of other Creatine Kinases offered on the market. For use, dissolve the lyophilized Creatine Kinase in MilliQ water at a concentration of 20 mg/mL. If you dissolve more of the enzyme than needed for your experiment, the solution should be dispensed into small aliquots and store at -80°C. Again, try to avoid repeated freezing and thawing of the Creatine Kinase solution. For our inhouse experiments, we commonly work only with fresh Creatine Kinase preparations.

Protocol

The WEPRO®8240 Series NMR Core Kits allow to perform protein expression reactions on different reaction scales. In this manual, we refer to a SMALL reaction scale of 240 µl (up to 100 reactions), a MEDIUM reaction scale of 1.2 ml (up to 20 reactions), and a LARGE reaction scale of 6 ml (up to 4 reactions). The reaction size is scalable and therefore other reaction sizes may also be used.

Note, the reagents provided with this kit can also be used to setup MEDIUM size reactions on a 1.2 ml scale (up to 16 reactions) and LARGE size reactions on a 6 ml scale (up to 3 reactions) on CellFree Sciences' fully automated Proteomist DTII protein synthesizer. Follow the manuals provided with the instrument for more details on how to use the reagents on the Proteomist DTII.

Bilayer translation reactions must be set up in flat-bottom vials! While individual flat-bottom vials are available, multiple reactions can be set up in multi-well plates. Multi-well plates are available with different dimensions to match the translation reaction size. It is preferable to have a larger surface between both layers to achieve higher protein yields.

Plate Size	Well Volume	Reaction Size	Name in Manual
96 wells	~ 360 µl	240 µl	SMALL
24 wells	~ 3.4 ml	1.2 ml	MEDIUM
6 wells	~ 16.8 ml	6 ml	LARGE


For first time use, we recommend including a positive control to make sure that the experiment has been set up correctly. Contact CFS on available control vectors that can be used in a separate expression experiment.

Successful protein expression should be confirmed before using any protein in your later experiments. Protein expression can be analyzed by SDS-PAGE to see whether a protein of the correct size has been made. It can be helpful to compare your protein in a crude expression reaction mixture to a negative control reaction prepared without added expression vector. The negative control reaction will only show the background proteins in the wheat germ extract. We advise to use known amounts of a BSA standard in the SDS-PAGE experiment to estimate protein yields. As an alternative, protein expression can also be confirmed by Western blotting using an antibody against the target protein or an affinity tag. When working with small protein amounts, Western blotting and labeling methods offer more sensitive protein detection than protein staining in SDS-PAGE gels. Those methods further provide background free detection as commonly proteins in the wheat germ extract should not be recognized by a specific antibody, nor is there any background when labeling proteins during expression experiments.

Time Requirements

Refer to the flowchart below on the estimated time per reaction step.

INSTRUCTION MANUAL

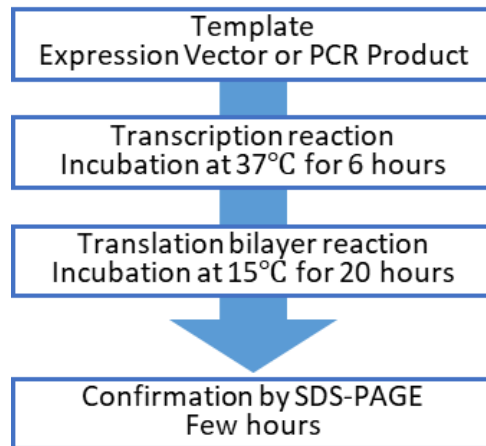

Transcription Reaction Using DNA Template

Per reaction perform the following steps to set up a transcription reaction on one of the reaction sizes:

1. Thaw your template DNA before the experiment. You need 1 µg (SMALL), 5 µg (MEDIUM) or 25 µg (LARGE) of purified plasmid DNA depending on the reaction scale (*1).
2. Take vials with 5xTranscription Buffer LM (yellow vial), NTP Mix (blue vial), RNase Inhibitor (orange vial) and SP6 RNA polymerase (green vial) out from storage at -80°C.
3. Thaw the reagents in a water bath at ~25°C. After thawing, spin the vials briefly to collect the entire volume at the bottom of the vial. Mix the reagents gently before use. Always keep reagents on ice.
4. Set up transcription experiment as shown in the table below for three different reaction sizes. Then mix gently by pipetting up and down.

Reagents	Volume SMALL	Volume MEDIUM	Volume LARGE	Final Concentration
Nuclease free water*	5.75 µl	28.8 µl	143.8 µl	-
5xTranscription Buffer LM	2 µl	10 µl	50 µl	1x
NTP Mix (25 mM)	1 µl	5 µl	25 µl	2.5 mM
RNase Inhibitor (80 U/µl)	0.125 µl	0.63 µl	3.13 µl	1 U/µl
SP6 RNA Polymerase (80 U/µl)	0.125 µl	0.63 µl	3.13 µl	1 U/µl
Plasmid (circular DNA, 1.0 µg/µl)	1 µl	5 µl	25 µl	100 ng/µl
Total	10 µl	50 µl	250 µl	

***Do not use DEPC treated water; we recommend using commercially available pure water.**

5. Incubate at 37°C for 6 hours in an incubator (*2).
6. After completion of the transcription reaction, leave reaction mixture at room temperature until later use in the translation reaction. Do not cool the reaction mixture, nor store it on ice.
7. Optionally, you can confirm the mRNA quality after the transcription reaction by agarose gel electrophoresis taking out 0.5 µl from the reaction mixture for loading onto the gel. Refer to a cloning handbook for more information on how to perform RNA gel electrophoresis. Successful RNA expression reactions yield in multiple transcripts that can be seen on the gel. Do not use an RNA preparation if you see an RNA smear because it indicates that your RNA had been degraded.

INSTRUCTION MANUAL

(Notes)

***1:** Commonly plasmid DNA prepared by a commercial DNA purification kit is suitable for use in protein expression experiments. Do not use plasmid DNA from alkaline lysis without further purification.

***2:** White precipitate may occur during incubation. This is magnesium pyrophosphate, which will not interfere with the following translation experiment. Use the whole reaction mixture including the precipitate in the next step.

Translation Reaction Using RNA Prepared from DNA Template

After completion of the transcription reaction, let the reaction mixture cool down to room temperature. Do not forcibly cool it down on ice or in a refrigerator. Resuspend the transcription mixture by pipetting gently up and down before use in translation reaction (*1).

Per reaction perform the following steps to set up translation reaction. Make sure to use information matching the transcription reaction size used in the previous step:

1. Take the vial with WEPRO®8240/WEPRO®8240G/WEPRO®8240H wheat germ extract (clear vial). Thaw the extract in a water bath at ~25°C. After thawing, briefly spin the vial with wheat germ extract to collect it at the bottom of the vial. Avoid excessive centrifugation of wheat germ extracts. Mix the extract gently before use. Keep wheat germ extract on ice at all time. After use, immediately freeze the wheat germ extract in liquid nitrogen and put it back into the freezer for storage at -80°C. The wheat germ extract loses its activity if not kept at -80°C!
2. Thaw Creatine Kinase (red vial) in a water bath at ~25°C. After thawing, briefly spin the vial with Creatine Kinase to collect it at the bottom of the vial. Mix the enzyme gently before use. We strongly recommend not to freeze/thaw Creatine Kinase as it will rapidly lose activity (*2).
3. Thaw 2xSUB-AMIX® SGC FOR NMR; refer to the table above on how to prepare the buffer.
4. Prepare “translation mixture” as indicated in the table below. Mix gently by pipetting up and down and avoid any bubbles.

Reagents	Volume SMALL	Volume MEDIUM	Volume LARGE	Final Concentration
Transcription Reaction	10 µl	50 µl	250 µl	1/2 volume
2xSUB-AMIX® SGC FOR NMR	19.2 µl	99.8 µl	499 µl	1x
Creatine Kinase (20mg/ml)	0.8 µl*	0.2 µl	1 µl	40 ng/µl
WEPRO®8240/WEPRO®8240G/WEPRO®8240H Extract	10 µl	50 µl	250 µl	120 OD/ml
Total	40 µl	200 µl	1000 µl	

***Dilute Creatine Kinase stock solution to 1 mg/ml with RNase free water when used for SMALL scale reaction setup!**

5. Dilute the 2xSUB-AMIX® SGC FOR NMR translation buffer with an equal volume of nuclease free water. Then add the required amount of 1xSUB-AMIX® SGC FOR NMR translation buffer to a flat-bottom vial or well. Refer to the table below on the required volumes for the different reaction sizes.

INSTRUCTION MANUAL

Reagents	Volume SMALL	Volume MEDIUM	Volume LARGE	Final Concentration
Translation Mixture	40 µl	200 µl	1000 µl	-
1xSUB-AMIX® SGC FOR NMR	200 µl	1.0 ml	5.0 ml	1x
Total	240 µl	1.2 ml	6 ml	

6. Carefully transfer the translation mixture from step 3 to the bottom of a single vial or well containing 1xSUB-AMIX® SGC FOR NMR translation buffer to form the bilayer reaction with the translation mixture in the lower layer and the 1xSUB-AMIX® SGC FOR NMR translation buffer in the upper layer. Refer to figure at the end of this section on how to setup a bilayer reaction: Go with the pipette tip to the bottom of the vial or well, and slowly release the translation mixture below the 1xSUB-AMIX® SGC FOR NMR translation buffer. Because of the yellowish color of the wheat germ extract, you can distinguish the translation mixture from the translation buffer. The wheat germ extract has a higher density than the buffer and therefore will form easily the lower layer.
- Do not mix the reagents in the vial or well by pipetting or any other means! It will reduce the yield of the reaction.**
7. Close the vial or seal the well to avoid evaporation.
8. Incubate at 15°C for 20 hours in an incubator. Be careful that the vial stably stands on a flat surface.
9. After completion of the translation reaction, mix the bilayer reaction gently by pipetting up and down.
10. Place reaction mixture on ice to protect your protein before later use.
11. Commonly some 2 to 3 µl of the crude translation reaction mixture are enough to detect the expressed protein on SDS-PAGE. It can be helpful to load different amounts of the crude reaction mixture to get a better estimation of the protein yields.

(Notes)

***1:** If you notice a white precipitate after the transcription reaction, resuspend the precipitate by pipetting gently up and down before mixing with the wheat germ extract. There is no need to remove the precipitate.

***2:** We recommend to always use fresh Creatine Kinase. Creatine Kinase is required for the energy supply of the translation reaction, and a loss of Creatine Kinase activity will reduce protein yields.

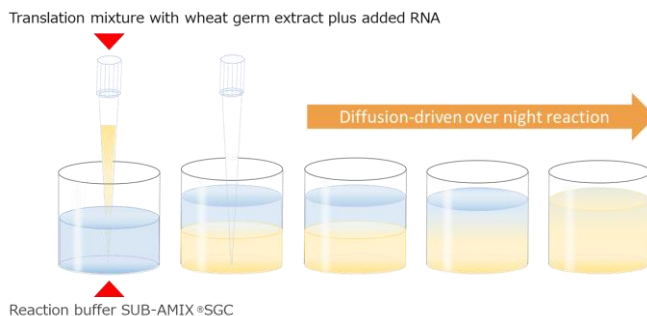


Illustration on how to setup bilayer reaction

Troubleshooting

The protein expression experiments require correct and accurate pipetting during reaction setup. Any mistake in the volumes added to the reactions, mixing the reagents, or forgetting any of the reagents will lead to wrong results. Therefore, carefully check the label for each reagent prior to starting the pipetting step.

- The experiment must be done under RNase-free conditions as any loss of the RNA template will prevent protein expression.
- Mark in your protocol each pipetting step you have completed.
- Change the pipetting tip after each pipetting step. Do not use the same pipetting tip to pipette different reagents or reaction mixtures. Always change the pipetting tip after use.
- Leaving out the plasmid template will always yield negative results. The same applies if there is a mistake in the expression vector, e.g. leaving out the starting ATG, forgetting a stop codon, or having a frame shift error.
- Confirm that your expression vector is correct and has a start and stop codon in line with the reading frame for the protein. Refer to our vector maps on more information on suitable sequencing primers to confirm the sequence of your expression vector.
- Confirm the DNA and RNA quality if the protein yields are low. Low RNA yields during the transcription reactions will also reduce the protein yields. Perform a phenol/chloroform extraction on the template DNA if RNA yields are low and ensure working under RNase-free conditions. An OD_{260/280} ratio of ~1.8 for your plasmid DNA preparations is commonly considered as pure enough for use in protein expression experiments. Lower ratios may indicate remaining proteins other contaminations absorbing near 280 nm. Note that the actual values for the OD_{260/280} ratio can vary from vector to vector as the actual OD values depend also on the nucleotide composition of your DNA vector. It is important to confirm the OD_{260/280} ratio for your vector DNA before use in protein expression experiments because low DNA purity prevents RNA and protein expression.
- We recommend to always using fresh Creatine Kinase. Do not freeze/thaw Creatine Kinase as it will rapidly lose activity. Creatine Kinase is required for the energy supply of the translation reactions, and a loss of Creatine Kinase activity will reduce protein yields.
- For the translation reaction, do not mix the two layers during setup of the bilayer translation reaction. Mixing both layers will sharply reduce the protein yields of a 20-hour translation reaction as the reaction will run dead within few hours. A slow mixing of both layers is required to maintain the translation reaction for up to 20 hours.
- It is possible to work with linear templates in cell-free protein expression experiments, which can be easily prepared by PCR methods. Working with PCR products, however, can reduce protein yields. PCR products are used to quickly find the best expression construct or to test expressing different protein fragments. If the yields obtained with PCR products are too low, it may be worthwhile to prepare an expression vector already for doing the test expression experiments. Always use expression vectors for large-scale protein production to have stable and reproducible conditions.
- If the protein is not expressed, check reaction conditions, reagents and DNA template in a small-scale expression reaction to confirm their integrity. If the results are unclear, you can check the

performance of the transcription and translation reactions separately to narrow down the problem. Use a positive control vector to make sure that all reagents work, and the experiment is done correctly. Reconsider the design of your expression template to improve protein yields if all the forgoing steps do not explain low protein yields.

- Make sure that the wheat germ extract was always kept frozen before use. Avoid repeated freeze/thawing; it will inactivate the extract.
- Store wheat germ extract at -80°C; storage at higher temperature will lead to low activity or even total loss of activity. Follow clearly the recommendations on reagent storage and handling.
- Keep all fractions during protein purification until you have confirmed the recovery of the purified proteins. If you are not able to recover the protein from the resin during the purification experiment, check whether the protein can be found in the flow through or the washing fractions.
- Some proteins may have special requirements and do not express well under standard conditions. Gather information on your target protein before the expression experiments to see whether additional considerations are needed. Contact us for more information on how to modify cell-free protein expression experiments.

Additional Information

Certain proteins may require changes to the expression reactions, where we have dedicated expression kits to prepare membrane proteins in the presence of added lipids or to prepare isotope-labeled proteins for use in MS or NMR studies. We can provide more information on the use of other additives such as detergents or ions in our cell-free protein expression system. Visit the homepage of CellFree Sciences for more information on other products and how to use our protein expression system.

Contact the technical support of CellFree Sciences for more information and further help. The contact information is given on the last page of the manual.

Customer Information

Label License Policy

Our products are covered by the US patents Nos. 7838640, 7981617, 8734856, and equivalent pending or granted patents in other countries. Please contact us at tech-sales@cfsciences.com if you intend to make commercial use of our expression system.

Commercial use includes:

- Using the product as a raw material in manufacturing process.
- Using the product for offering paid services.
- Reselling the product or any of its components.

All our products are for RESEARCH USE ONLY. NOT FOR HUMAN OR ANIMAL THERAPEUTIC OR DIAGNOSTIC USE.

Trademarks

ENDEXT[®], WEPRO[®] and SUB-AMIX[®] are registered trademarks of CellFree Sciences Co., Ltd.

Others

All product specifications and information in the manual may be changed without prior notice.

Contact

CellFree Sciences Co., Ltd.

Yokohama Head Office

Yokohama Bio Industry Center, 1-6 Suehiro-cho, Tsurumi-ku, Yokohama, 230-0045 JAPAN

TEL : +81-45-345-2625 FAX : +81-45-345-2626

E-mail: tech-sales@cfsciences.com

Follow us on Twitter (@CFSciences) and LinkedIn for up-to-date information on new publications, products, and the use of our expression system.

CellFree Sciences is certified under ISO 9001:2015. - Our products and services are for research purpose only. - Refer to our homepage at <http://www.cfsciences.com/eg/> for more information on our products and services, or contact us directly at tech-sales@cfsciences.com.

Appendix

SUB-AMIX® SGC FOR NMR (¹⁵N)

List of labeled amino acids provided with the ¹⁵N versions of the WEPRO®8240 Series NMR Core Kit:

Amino Acid	Label	Molecular Weight
L-ASPARTIC ACID	15N, 98%	134.1
L-ARGININE:HCL	15N4, 98%	214.64
L-CYSTEINE	15N, 98%	122.15
L-GLUTAMIC ACID	15N, 98%	148.12
L-TRYPTOPHAN	15N2, 98%	206.21
L-TYROSINE	15N, 98%	182.18
L-ALANINE	15N, 98%	90.09
L-ASPARAGINE:H2O	15N2, 98%	152.12
L-GLUTAMINE	15N2, 98%	147.13
GLYCINE	15N, 98%	76.06
L-HISTIDINE:HCL:H2O	15N3, 98%	212.61
L-ISOLEUCINE	15N, 98%	132.17
L-LEUCINE	15N, 98%	132.17
L-LYSINE:2HCL	15N2, 98%	221.09
L-METHIONINE	15N, 98%	150.21
L-PHENYLALANINE	15N, 98%	166.18
L-PROLINE	15N, 98%	116.12
L-SERINE	15N, 98%	106.09
L-THREONINE	15N, 98%	120.11
L-VALINE	15N, 98%	118.14

INSTRUCTION MANUAL

SUB-AMIX® SGC FOR NMR (¹³C/¹⁵N)

List of labeled amino acids provided with the ¹³C/¹⁵N versions of the WEPRO® 8240 Series NMR Core Kit:

Amino Acid	Label	Molecular Weight
L-ASPARTIC ACID	U-13C4, 97-99%; 15N, 97-99%	138.07
L-ARGININE:HCL	U-13C6, 97-99%; 15N4, 97-99%	220.59
L-CYSTEINE	U-13C3, 97-99%; 15N, 97-99%	125.13
L-GLUTAMIC ACID	U-13C5, 97-99%; 15N, 97-99%	153.09
L-TRYPTOPHAN	U-13C11, 97-99%; 15N2, 97-99%	217.13
L-TYROSINE	U-13C9, 97-99%; 15N, 97-99%	191.12
L-ALANINE	U-13C3, 97-99%; 15N, 97-99%	93.07
L-ASPARAGINE:H2O	U-13C4, 97-99%; 15N2, 97-99%	156.09
L-GLUTAMINE	U-13C5, 97-99%; 15N2, 97-99%	153.1
GLYCINE	U-13C2, 97-99%; 15N, 97-99%	78.05
L-HISTIDINE:HCL:H2O (<5% D)	U-13C6, 97-99%; 15N3, 97-99%	218.57
L-ISOLEUCINE	U-13C6, 97-99%; 15N, 97-99%	138.12
L-LEUCINE	U-13C6, 97-99%; 15N, 97-99%	138.12
L-LYSINE:2HCL	U-13C6, 97-99%; 15N2, 97-99%	227.05
L-METHIONINE	U-13C5, 97-99%; 15N, 97-99%	155.17
L-PHENYLALANINE	U-13C9, 97-99%; 15N, 97-99%	175.12
L-PROLINE	U-13C5, 97-99%; 15N, 97-99%	121.09
L-SERINE	U-13C3, 97-99%; 15N, 97-99%	109.06
L-THREONINE	U-13C4, 97-99%; 15N, 97-99%	124.08
L-VALINE	U-13C5, 97-99%; 15N, 97-99%	123.1